

### Fixatives

GLUTARALDEHYDE	
2.5% glutaraldehyde/0.1 M buffer	5 ml 25% glutaraldehyde 20 ml distilled water 25 ml 0.2M buffer
2.5% glutaraldehyde/0.1 M buffer + 1% tannic acid	0.5 g tannic acid dissolved in 20 ml distilled water 5 ml 25% glutaraldehyde 25 ml 0.2M buffer
OSMIUM TETROXIDE	
1% OsO <sub>4</sub> / 0.1 M buffer	5 ml 4% OsO <sub>4</sub> (EMS-19150) 5 ml distilled water 10 ml 0.2 M buffer